

Unrivalled RNA Purity

TRIsure™

- Ready-to-use solution for Isolation of premium quality Total RNA
- Column-free Total RNA Isolation, optimized for maximum RNA yield
- Suitable for a wide variety of cells and tissues including bacteria and plants

TRIsure™ is a ready-to-use reagent for the isolation of high quality total RNA from diverse biological materials, including animal tissues, cultured cells, bacterial cells, as well as plant tissues rich in polysaccharides and proteoglycans. TRIsure maintains the integrity of the extracted RNA (high RIN and 28S/18S ratio) (Fig. 1), while disrupting cells and dissolving cell components. The reagent combines a proprietary blend of phenol and other components for optimal results.

The biological sample is homogenized or lysed in TRIsure and then separated into organic and aqueous phases. The RNA remains in the aqueous phase and is subsequently recovered by precipitation with isopropyl alcohol. High yield (Table 1) and high quality RNA, with virtually no genomic DNA contamination, is extracted (Fig. 2). A volume of 1ml of TRIsure is sufficient to isolate total RNA from 1×10^7 cells or 100mg of tissue. The isolation method is rapid and straightforward (Fig. 3).

Using intact RNA as provided by TRIsure, is a key element in obtaining reliable gene expression data in downstream applications, such as cDNA synthesis, real-time RT-PCR, microarray, hybridization assays and *in vitro* translation (see product citations).

Table 1. Expected yield of RNA from various samples using TRIsure

Sample type	Sample quantity	Expected yield
Cultured epithelial cells	1×10^6	8-15µg
Cultured fibroblasts	1×10^6	20-25µg
Mouse kidney tissue	1mg	2-5µg
Mouse liver tissue	1mg	5-10µg

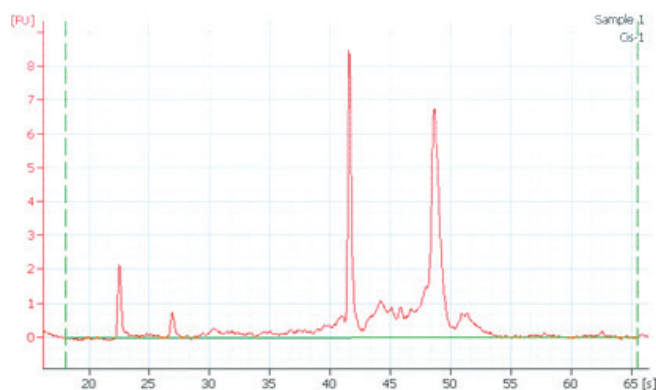


Fig. 1 Isolation of high quality RNA.

A log phase culture of *Bacillus subtilis* was pre-treated with Bacterial Enhancement Reagent, followed by isolation of RNA using TRIsure. The RNA was analyzed using Bioanalyzer 2100 (Agilent Technologies) and was found to be of high quality and purity.



United Kingdom
Tel: +44 (0)20 8830 5300

Germany
Tel: +49 (0)3371 681 229

USA
Toll Free: 888 257 5155

Australia
Tel: +61 (0)2 9209 4180

www.bioline.com/rna



A Meridian Life Science® Company

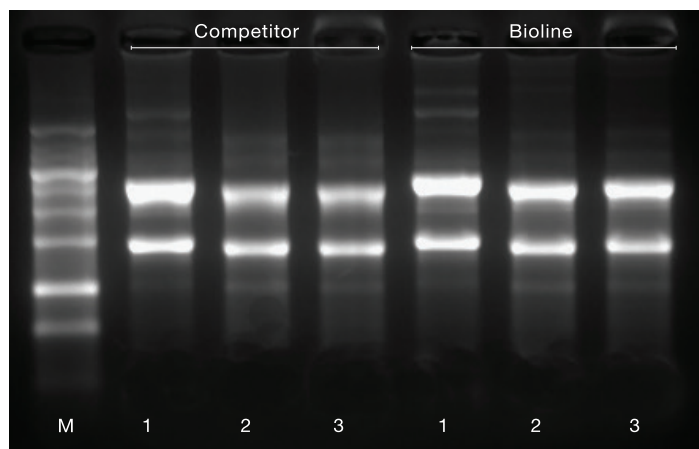


Fig. 2 High quality and yield of RNA extracted using TRI[®]sure.

RNA extracted from 3T6 cells and mouse tissue, using TRI[®]sure and Competitor reagent.

Lane 1: 4µg of total RNA from 3T6 cells

Lane 2: 4µg of total RNA from mouse kidney tissue

Lane 3: 4µg of total RNA from mouse liver tissue


M: RNA Ladder

Product Citations

1. Chincinska, I.A., *et al. Plant Physiol.* **146**,515-528 (2008).
2. Hedtke, B., *et al. NAR* **37**(11), 3739-3746 (2009).
3. Tsuda, M., *et al. J. Cell. Mol. Med.* **13**, 412-603 (2009).
4. Schroeckh, V., *et al. PNAS* **106**, 14558-14563 (2009).
5. Yamazaki, S., *et al. Biochim. Biophys. Acta* **1779**, 2108-2114 (2009).
6. Gollan, P. J., *et al. Physiologia Plantarum* **143**(4), 385-395 (2011).
7. Wanger, A. E., *et al. BMC Comp. Alt. Medicine* **11**, 1 (2011).
8. Yoshida, A., *et al. Plant J.* **70**(2), 327-339 (2012).



Fig. 3 TRI[®]sure Protocol.

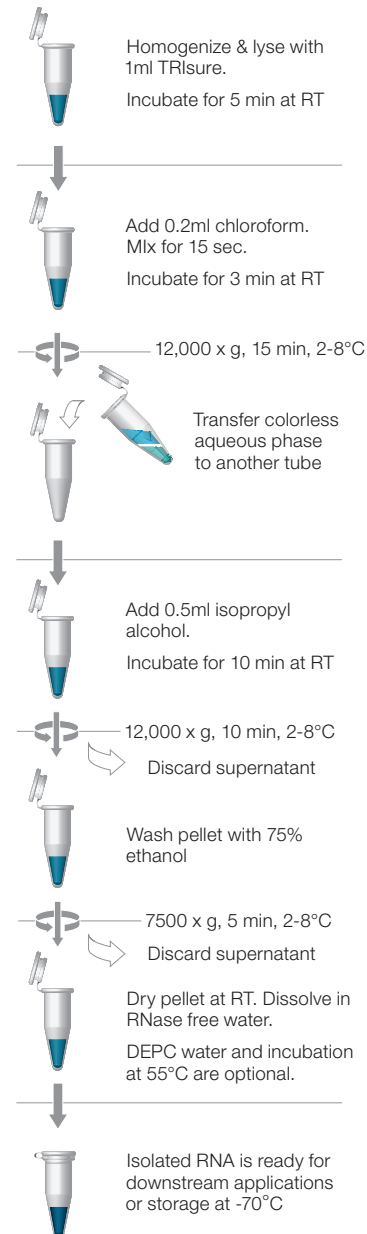

Ordering Information

PRODUCT	PACK SIZE	CAT NO.
TRI [®] sure	100ml	BIO-38032
TRI [®] sure	200ml	BIO-38033

PS0612V3



United Kingdom
Tel: +44 (0)20 8830 5300

Germany
Tel: +49 (0)3371 681 229

USA
Toll Free: 888 257 5155

Australia
Tel: +61 (0)2 9209 4180

www.bioline.com/rna



A Meridian Life Science® Company